

Ethanolamines/Glycol Ethers

Tier 3 Procedure

Section Number: 606
Issue Number: 1.5
Original Date: 12-17-91 **Author:** Paula Bourgeois
Revision Date: 03-11-98
Title: Preparing T-240 Overheads Sample: Gas Chromatograph #'s 7 & 9
References: Glossary Of Equipment Identification (Text # 114)

=====

606. Preparing T-240 Overheads Sample For Gas Chromatograph #'s 7 And 9

1.0 Objective:

1.1 To ensure sample is prepared consistently from person to person for more accurate analysis.

2.0 Responsibility:

2.1 It is the responsibility of each operator performing this task to comply with this procedure.

3.0 Procedure:

3.1 Draw up 1.0cc of sample using a 1.0cc plastic disposable syringe.

3.2 Inject the sample into a GC vial.

3.3 Cap the vial with an aluminum seal cap and the crimper.

3.4 Set up the ethanolamines hp 5890 GC using the instructions on the multichron system.

4.0 Revisions:

Revision No. 1	02-19-93	Yearly Review
Revision No. 2	12-21-94	Yearly Review
Revision No. 3	02-24-95	January/Yearly review
Revision No. 4	12-29-95	Yearly Review
Revision No. 5	03-11-98	Modified 3.4