

CONFIDENTIAL:
Subject to Protective Order
of 14th Judicial District Court
No. 91-1145

PLANT LAB EQUIPMENT MAINTENANCE

INTEGRATORS:

WEEKLY----- VACUUM FILTERS, LOCATED BELOW KEYBOARD
ON FRONT OF 4100'S AND 4200'S.
----- NEVER ALLOW ANY SAMPLES WITH A CORROSIVE
NATURE NEAR THE INSTRUMENT.

G.C.'S:

WEEKLY----- VACUUM AIR INTAKE VENTS.

EVERY

SHIFT-----

CHANGE SEPTUM.

----- CHECK GAS BOTTLES, CHANGE WHEN BELOW
500 psi.

--- CHECK OPERATION OF VACUUM PUMP USED TO
CLEAN SYRINGES.

--- NEVER ALLOW ANY SAMPLE WITH A CORROSIVE
NATURE NEAR THE INSTRUMENT.

WET TEST INSTRUMENTS:

DAILY----- WIPE SURFACE WITH A CLEAN, DAMP TOWEL.

----- CHECK REFERENCE SOLUTION IN pH
ELECTRODE.

--- NEVER ALLOW ANY SAMPLE WITH A CORROSIVE
NATURE NEAR THE INSTRUMENT.

SL 002190